



A SIGN OF QUALITY

Set Up Instructions for Vaccine Refrigerators

Dear Customer,

Thank you for choosing our refrigerator, to ensure that cabinet is installed correctly please follow the following steps.

- a. Unpack the refrigerator and place in position
- b. Ensure that refrigerator is level, if not you may adjust front levelling feet only
- c. Ensure there is 60mm ventilation space at rear and 30mm on sides of refrigerator for proper air flow.
- d. Position shelves at the desired levels that you require.
- e. Plug refrigerator into your 240 volts outlet and let run for 2 hours during the period the refrigerator would have cycled a number of times. You are now ready to place your products on shelves.
- f. Cabinet temperature will vary from when the motor stops and starts. This is only air temperature and not product temperature.
- g. Temperature reading should only be done as follows, reset controller at end of day and do not open door again, next morning take your high and low read out before you open door.

Preventative Maintenance

It is important that the condenser on the back of the cabinet is cleaned every 3 months, simply by brushing the condenser in a vertical motion with a small brush or vacuum cleaner.

However this will also depend on how dusty the area is. You may be able to do it every 6 months.

Refrigeration is fan forced to ensure even temperature throughout the cabinet with auto defrost and water disposal.